UPGRADE KIT RTD-502 / CK-7

Instructions for installing the Upgrade Kit "RTD-502 / CK-7"

Upgrade Kit contains:

1. 1 electronics insert RTD-502, interchangeable
2. 1 new resistive electrolyte measuring cell CK-7 with "Novalyte" technology
3. 1 sticker red "RTD-502/CK-7" for back panel placing

This upgrade-kit can be installed at a customer.

Upgrade Kit 230V : Art. No. **111 8687**.
 Upgrade Kit 100/115V : Art. No. **111 8688**

Installation procedure on site

by customer or by our representative's service department

Step 1, electronics insert RTD:

- Disconnect instrument from mains.
- Remove the 2 screws at top and bottom of the front panel of the electronics RTD-500.
- Pull out the RTD-500 slowly.
- Carefully completely push in the new RTD-502. Then lock it with the 2 screws.

Step 2, measuring head BSK:

- Remove the measuring cell CK from the head BSK. This is quite simple, when using the Novasina tweezers specially developed for that purpose: optional art. 1111218. Now, following the instructions in the operating manual of your instrument, very carefully insert the new cell CK-7 (flag towards white dot).

Step 3, sticker:

- On the back panel, place the red sticker "RTD-502/CK-7" near the sensor cable connector. This is to show, that the instrument now is equipped with a cell type CK-7.

Step 4, calibration:

- As after every cell replacement, a new calibration is necessary. A factory precalibration at 6 points from 11 to 98%rh has already been done!

!! After modification: return the RTD-500 for crediting !!

Enclosure to the operation manual

UPGRADE KIT RTD-502 / CK-7

New
RTD-502 (Software V3.13 or higher) / CK-7

 measuring range: **0.03 ... 1.00aw** (3 ... 100%rh).

Operating temperature range: 0°C ... 50°C.

The display will show "low hum", if the measurement goes below ca. 0.025aw.

Modification of the calibration procedure compared to the RTD-500-electronics:

SW versions V2.0-2.20:

- table with all calibration values: if a point is not calibrated, the display now shows " -.- ", because an internal reference value of " 0.0 " is possible at 4%.
- the text "low cal" was removed.
- **New:** automatic calibration points at 6% and at 4%rh.

SW version V3.12:

- **New:** additional automatic calibration point at 4%rh.

Calibration of an instrument, following the operation manual:

- the first = basic calibration point is still with SAL-T 75%. At this point, also the cell characteristics ("HC") is determined.
- the sequence of the following points can freely be chosen.
- special procedure at 11%: when placing a Novasina SAL-T 11%, a SAL-T 6%, or a 4% standard (CsF made by the user, not available from Novasina), and trying to calibrate, the menu is displaying the following question: "cal 11?". This question must be answered by confirming (if salt 11 is in the chamber), or going with up/down button to "cal 6" or "cal 4", depending on the humidity standard in the chamber. Then confirm.

The "calibration protocol" was also modified to show all 8 calibration points.

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